

## Introduction

The STEMPRO® Osteogenesis Differentiation Kit has been developed for the osteogenic differentiation of mesenchymal stem cells (MSCs) in tissue-culture vessels. The kit contains all reagents required for inducing MSCs to be committed to the osteogenesis pathway and generate osteocytes. Using STEMPRO® Osteogenesis Differentiation Kit in combination with STEMPRO® MSC SFM or MesenPRO RS™ Medium provides a standardized culture workflow solution for MSC isolation, expansion, and differentiation into matrix-forming osteocytes.

Description	Cat. no.	Size	Storage	Shelf Life
STEMPRO® Osteogenesis Differentiation Kit	A10072-01	1 kit		—
Contains:				
STEMPRO® Osteocyte/Chondrocyte Differentiation Basal Medium	A10069-01	100 mL	2 to 8°C (protect from light)	12 months
STEMPRO® Osteogenesis Supplement	A10066-01	10 mL	-5 to -20°C (in the dark)	12 months


## Intended Use

For research use only (RUO). **Caution:** Not intended for human or animal diagnostic or therapeutic uses.


## Characteristics

The STEMPRO® Osteogenesis Differentiation Kit has been extensively tested and proven to have the following characteristics:

- Contains all components required to reliably and reproducibly induce MSCs into the osteogenic lineage.
- Demonstrated to robustly induce osteogenesis in adipose tissue-derived stem cells (STEMPRO® Human Adipose-Derived Stem Cell Kit, Cat. nos. R7788-110 and R7788-115).
- Classical staining methods demonstrate differentiation of MSCs into osteoblasts and osteocytes (Figure 1).
- Immunocytochemistry methods demonstrate expression of relevant osteocyte biomarkers (Figure 2).

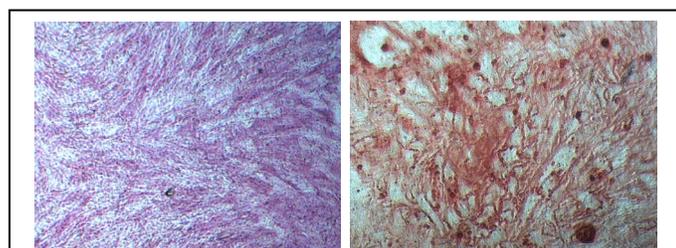

**Figure 1:** Analysis of MSCs cultured in STEMPRO® Osteogenesis Differentiation Medium demonstrated differentiation into osteogenic lineages by alkaline phosphatase staining and Alizarin Red S staining.

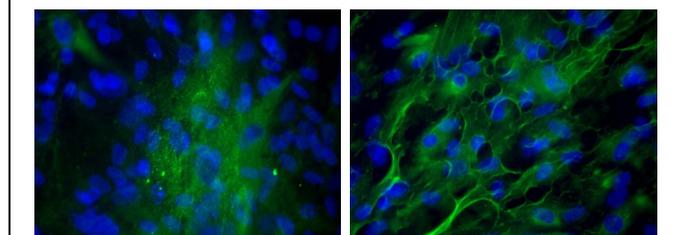

**Figure 2:** Analysis of MSCs cultured in STEMPRO® Osteogenesis Differentiation Medium demonstrated differentiation into osteogenic lineages by bone sialoprotein and osteopontin immunostaining.

## Storage and Handling

- STEMPRO® Osteogenesis Supplement is supplied frozen. Thaw prior to use, as described in **Media Preparation**, next page.
- Thawed STEMPRO® Osteogenesis Supplement is stable up to at least one month at 2 to 8°C. Supplement can be refrozen in desired volumes and stored at -5°C to -20°C. **Avoid multiple freeze thaw cycles** of supplement.
- Store prepared Complete STEMPRO® Osteogenesis Differentiation Medium at 2 to 8°C in the dark. Complete medium is stable up to at least one month at 2 to 8°C.

## Important Guidelines for Osteogenesis Differentiation

To obtain optimal osteogenic differentiation with STEMPRO® Osteogenesis Differentiation Medium, follow these guidelines:

- **Expansion culture:** Primary MSC isolates should be expanded with STEMPRO® MSC SFM or MesenPRO RS™ Medium in T-75 or T-225 flasks. Standard growth media of DMEM+10% MSC Qualified FBS has been successfully tested. We recommend refeeding the cultures every 2 to 3 days and passaging every 5 to 7 days.
- **Passaging:** We strongly recommend using low-passage MSCs (<8 to 10 passages). Continuously passaged MSCs will gradually lose their multipotency with increased passage number (>10 passages).
- **Harvesting:** We recommend using TrypLE™ Express for enzymatically treating and harvesting MSCs. TrypLE™ Express is a recombinant protease that has been demonstrated to be gentle on MSCs. Overexposure to trypsin will lead to reduced MSC viability and expansion.
- **Timing of passaging:** It is critical to not let passaged MSCs become completely confluent, as it can reduce multipotency of MSCs. Passage cultures when they reach 60 to 80% confluency, cell viability is at least 90%, and the growth rate is in mid-logarithmic phase.
- **Seeding density:** For expansion, we recommend a seeding density of  $3 \times 10^3$  to  $5 \times 10^3$  viable cells/cm<sup>2</sup> with MesenPRO RS™ Medium or  $1 \times 10^4$  viable cells/cm<sup>2</sup> with STEMPRO® MSC SFM.
- **Confluency:** Expanding MSCs in growth medium for 2 to 4 days before refeeding with Complete Osteogenesis Differentiation Medium can enhance osteogenesis.

## Certificate of Analysis

The Certificate of Analysis (CofA) provides quality control information for this product. The CofA is available on our website at [www.invitrogen.com/cofa](http://www.invitrogen.com/cofa), and is searchable by product lot number, which is printed on the box.

## Physical Conditions for Osteogenesis Culture

**Media:** STEMPRO® Osteogenesis Differentiation Medium

**Cell Line:** Human mesenchymal stem cells

**Incubator:** 36 to 38°C, humidified atmosphere of 4 to 6 % CO<sub>2</sub> in air

**Culture Conditions:** Adherent; ensure proper gas exchange and minimize exposure to light

**Recommended Culture Vessels:** 12-well tissue-culture plates, 16-well CultureWell™ slides, 96-well tissue-culture plates, or 75-cm<sup>2</sup> tissue-culture flasks

## Media Preparation

**Complete Osteogenesis Differentiation Medium:** Thaw supplement at 4°C, room temperature, or in a 37°C water bath, and prepare as below. Store complete medium at 2 to 8°C in the dark.

Osteogenesis Differentiation Medium	Final Conc.	For 100 mL
STEMPRO® Osteocyte/Chondrocyte Differentiation Basal Medium	1X	90 mL
STEMPRO® Osteogenesis Supplement	1X	10 mL
Gentamicin (10 mg/mL)	5 µg/mL	50 µl

**MSC Growth Medium:** Prepare as below.

MSC Growth Medium	Final Conc.	For 500 mL
DMEM low glucose		445 mL
MSC-qualified FBS	10%	50 mL
GLUTAMAX™ -I (200 mM)	2 mM	5 mL
Gentamicin (10 mg/mL)	5 µg/mL	250 µl

## Osteogenesis Differentiation

1. Observe cell monolayer from basal cultures expanded in STEMPRO® MSC SFM, MesenPRO™ RS medium, or standard growth medium (DMEM+10% FBS) to ensure mid-log growth phase confluence (60 to 80%). Aspirate medium and floating cells from culture flask and discard.
2. Add 5 to 10 mL DPBS. Gently rinse cell monolayer.
3. Remove DPBS, add 5 to 7 mL of pre-warmed TrypLE™ Express to flask and completely coat culture surface. Incubate for 5 to 8 minutes at 36 to 38°C or until cells have fully detached.
4. Gently pipet detached cells into a single cell solution and verify on inverted microscope.
5. Remove cell suspension from flask, transfer into a centrifuge tube, and pellet cells at 100 x g for 5 to 10 minutes.
6. Determine cell viability and total cell density using Trypan Blue Stain and electronic (*i.e.*, Coulter Counter) or manual (*i.e.*, hemocytometer) cell counting method.
7. Resuspend pellet in an appropriate volume of pre-warmed MSC Growth Medium (see **Media Preparation**).
8. Seed MSCs into culture vessels at 5 x 10<sup>3</sup> cells/cm<sup>2</sup>. For classical stain differentiation assays, seed into a 12-well plate. For gene expression profile studies, seed into a T-75 flask. For immunocytochemistry studies, seed into a 16-well CultureWell™ chambered coverglass or 96-well plate.
9. Incubate in MSC Growth Medium at 36 to 38°C in a humidified atmosphere of 4 to 6% CO<sub>2</sub> for a minimum of 2 hours up to 4 days.
10. Replace media with pre-warmed Complete Osteogenesis Differentiation Medium and continue incubation. MSCs will continue to expand as they differentiate under osteogenic conditions. Refeed cultures every 3 to 4 days.
11. After specific periods of cultivation, osteogenic cultures can be processed for alkaline phosphatase staining (7 to

14 days) or Alizarin Red S staining (>21 days; see below for method), gene expression analysis, or protein detection.

## Alizarin Red S Stain Analysis

1. After 21 days or longer under differentiating condition, remove media from 12-well plate and rinse once with DPBS. Fix cells with 4% formaldehyde solution for 30 minutes.
2. After fixation, rinse wells twice with distilled water and stain cells with 2% Alizarin Red S solution (pH 4.2) for 2 to 3 minutes.
3. Rinse wells three times with distilled water, visualize under light microscope and capture images for qualitative or quantitative analysis.


## Additional Products

Some products are recommended but not supplied in the kit. See below for ordering information.

Product	Size	Cat. no.
STEMPRO® MSC SFM	500 mL	A10332-01
STEMPRO® Human Adipose-Derived Stem Cell Kit	1 kit	R7788
STEMPRO® Adipogenesis Differentiation Kit	1 kit	A10070-01
MesenPRO RS™ Medium	500 mL	12746
FBS, MSC-Qualified (non-US)	100 mL	12662
GLUTAMAX™ -I	100 mL	35050
Gentamicin (10 mg/mL)	10 mL	15710
TrypLE™ Express	100 mL	12604
DPBS without Ca <sup>++</sup> and Mg <sup>++</sup>	500 mL	14190
Mouse anti-Osteocalcin	100 µg	33-5700
Rabbit anti-Osteopontin	100 µg	42-7701
CultureWell™ chambered coverglass	1 pack	C-37005
Trypan Blue Stain	100 mL	15250

## Explanation of Symbols and Warnings

The symbols present on the product label are explained below:

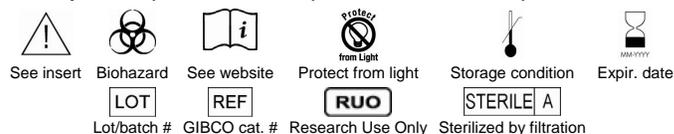

## Purchaser Notification

This product is covered by Limited Use Label Licenses (see our web-site [www.invitrogen.com](http://www.invitrogen.com), or the Invitrogen catalog). By the use of this product you accept the terms and conditions of all applicable Limited Use Label Licenses.

**Limited Use Label License No. 5: Invitrogen Technology:** The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. For products that are subject to multiple limited use label licenses, the most restrictive terms apply. Invitrogen Corporation will not assert a claim against the buyer of infringement of patents owned or controlled by Invitrogen Corporation which cover this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Licensing Department, Invitrogen Corporation, 1600 Faraday Avenue, Carlsbad, California 92008. Phone (760) 603-7200. Fax (760) 602-6500. Email: [outlicensing@invitrogen.com](mailto:outlicensing@invitrogen.com).

GIBCO® and STEMPRO® are registered trademarks and MesenPRO RS™, TrypLE™, and GLUTAMAX™ are trademarks of Invitrogen Corporation. Culturewell™ is a trademark of Grace Bio-Labs, Inc.

## Technical Support

Worldwide email: [techsupport@invitrogen.com](mailto:techsupport@invitrogen.com). Toll-free U.S. phone support: 1 800 955 6288. For additional country-specific support, visit our website at [www.invitrogen.com/contacts](http://www.invitrogen.com/contacts).